

Enzymax LLC, 870 Corporate Dr., Suite 201; Web: www.enzymax.net; Tel: 859-219-8482; Fax: 859-219-0653

EZshredder for simple and quick homogenization of cell and tissue lysates

Cat. #	Description	Quantity
EZC109S	EZshredder for Tissue & Cell-lysate homogenization (replace syringe and needle) in nucleic acid minipreps and protein purification to remove insoluble debris and reduce viscosity (Equivalent to QIAshredder)	50
EZC109L	EZshredder for Tissue & Cell-lysate homogenization (replace syringe and needle) in nucleic acid minipreps and protein purification to remove insoluble debris and reduce viscosity (Equivalent to QIAshredder)	250

All Bulk columns come with 2ml collection tubes with caps

Description:

EZshredder column provides an easy and efficient spin column method for homogenization of cell or tissue lysates. The column contains a biopolymer-shredding membrane which is in a microcentrifuge spin-column format. Here are the Highlights for the column:

Highlights/Applications:

- □ Quick and easy spin to replace syringe and needle homogenization
- □ Prevent sample material loss by simply loading the sample to column and spin to collect the flow-through
- \Box No cross-contamination between samples unless using the same column for different samples \Box Spin to remove the insoluble debris
- □ Reduce viscosity of samples for downstream RNA or miRNA preparation (with different nucleotide biding columns) or protein purification to obtain high quality and high yield products □ Shear genomic DNA (<20kb) to reduce viscosity for protein sample loading on PAGE gel
- □ Shear genomic DNA (<20kb) to reduce viscosity for protein sample loading on PAGE gel □ Filter Disc does not bind to DNA/RNA or Proteins
- THE DISC GOES HOLDING TO DIVA/KINA



EZshredder Spin Column (for Tissue & Cell- lysate homogenization)

How to use it:

- Load cell or tissue lysate into the EZshredder column with a collection tube.
- Spin at full speed in a microcentrifuge.
- Collect the flow-through in the tube.
- Samples are ready for downstream application, eg, RNA or miRNA mini/midi preparation or protein purification.